asecos®

SP  
LINE

### USER MANUAL

Safety storage cabinets according to  
SP Method 2369 version 6 (2018)

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#### 5 YEARS WARRANTY\*

\*Upon conclusion of an asecos service and maintenance agreement (**BasicPlus tariff**) with a fixed term of 5 years, you will get a warranty extension for a maximum of 5 years for your safety storage cabinet.

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For detailed information about our warranties please visit  
[www.asecos.com/service-conditions](http://www.asecos.com/service-conditions)

# SP LINE

SP30.145.085.NE.WDFW  
SP30.145.085.NE.WDFWR

***Models prepared for  
technical ventilation***

SP30.145.085.EA.WDFW  
SP30.145.085.EA.WDFWR

SP30.195.055.NE.WDFW  
SP30.195.055.NE.WDFWR

***Models prepared for  
technical ventilation***

SP30.195.055.EA.WDFW  
SP30.195.055.EA.WDFWR

SP30.195.085.NE.WDFW  
SP30.195.085.NE.WDFWR

***Models prepared for  
technical ventilation***

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asecos GmbH

## **YOUR PERSONAL DOCUMENTATION TO THE asecos SAFETY CABINET**

Dear Customer,

you have made a decisive investment in safety for your company by purchasing this asecos safety storage cabinet. You now own an innovative product made of top-quality materials guaranteeing the highest quality standards.

asecos safety storage cabinets have complete authorisation documents. We archive the authorisation documents for every individual cabinet, keeping them ready for you should you ever need them (e.g. for a works inspection or similar). Simply request them using this form.

Tear off or copy that page and return to us by fax with your address and serial number of the cabinet on it.

Yours sincerely

asecos GmbH

<b>Firma</b> • company	
<b>Abteilung</b> • department	
<b>Name</b> • name	
<b>Straße</b> • street	
<b>PLZ</b> • postal code	
<b>Ort</b> • city	
<b>Schrank</b> • cabinet <b>Seriennummer(n)</b> • serial number(s)	

# USER MANUAL

Dear customer,

Thank you very much for purchasing a safety cabinet from our company, with which you have made a decisive investment for safety in your home. With our safety cabinets the storage of hazardous substances becomes safe and comfortable for you.

Please read these operating instructions very carefully. Get to know the advantages and easy handling of our safety cabinets in detail. This will simplify your daily handling of hazardous substances

Thanks a lot  
Your asecos team

---

## USER MANUAL



<b>1. NOTES • GUIDELINES • GUARANTEE</b> .....	<b>5</b>
1.1. General safety notes .....	5
1.2. Guarantee .....	5
1.3. Cabinet details .....	5
<b>2. TRANSPORT</b> .....	<b>6</b>
2.1. Tilting the cabinet .....	6
2.2. In-plant transport .....	6
<b>3. ERECTION • COMMISSIONING • FUNCTION</b> .....	<b>6</b>
3.1. Alignment of the cabinets .....	6
3.2. Commissioning .....	7
3.3. Interior lighting (optional) .....	7
<b>4. CLOSING MECHANISM</b> .....	<b>7</b>
<b>5. INTERIOR EQUIPMENT</b> .....	<b>8</b>
5.1. Bottom collecting sump .....	8
5.2. Shelves (height-adjustable) .....	8
<b>6. STORAGE</b> .....	<b>9</b>
<b>8. FIRE • DISPOSAL</b> .....	<b>9</b>
8.1. Opening the cabinet after the fire .....	9
8.2. Disposal .....	10
<b>9. SAFETY CHECKS</b> .....	<b>10</b>
9.1. General information .....	10
9.2. Contact .....	10
<b>10. TECHNICAL DATA</b> .....	<b>11</b>
<b>11. TECHNICAL DRAWING</b> .....	<b>11</b>
<b>12. CABINET MARKING</b> .....	<b>13</b>

## 1.1. GENERAL SAFETY NOTES

- Observe applicable statutes and regulations, and the notes in these operating instructions, when handling hazardous materials.
- Follow the instructions in SP Method 2369 for the storage of flammable products in sales rooms
- Work on the electrical system is to be carried out only with the power turned off and only by qualified electricians – refer here to the regulations of the local electricity supply company.
- The on-site installation conditions are to be observed
- The instructions of the supervisory engineering department must be followed.
- Observe accident prevention regulations and workplace ordinance.
- Ensure that the necessary safety checks are only carried out by authorised staff using original spare parts.
- The pivoting range of the doors must always be kept clear, doors must be kept closed
- Observe the upper limits for stored quantities, loading etc.
- Containers with a volume larger than that of the spill containment sump may not be placed inside the cabinet; spilt hazardous materials are to be collected immediately and removed.
- Containers with aggressive chemicals (acids/alkalis) are to be kept in special cabinets for acids and alkalis or on shelves for acids and alkalis as well as in cabinets with metal-free interior equipment.
- Before storing such materials, check that the cabinet's surface is resistant to the chemicals.
- Before the initial commissioning, the safety storage cabinet is to be examined by the user for possible damage (see point 3.2)

### Set-up requirements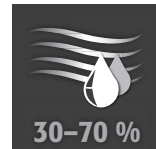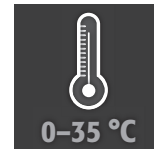

## 1.2. GUARANTEE

The guarantee for this product is agreed between you (the customer) and your dealer (the seller). As the manufacturer, asecos guarantees the products listed in the operating instructions for a period of 24 months from the date of delivery.

All model safety equipment are subject to a compulsory annual inspection by specialised staff authorised by the manufacturer. Otherwise the customer's guarantee claim against the manufacturer expires.

## 1.3. CABINET DETAILS

**You will find a complete model overview at the beginning of the operating instructions**

Cabinet data	⇒ logbook (included with the cabinet)
Technical drawing	⇒ page 11
Technical data	⇒ page 11

### SP-LINE

These models are tested, certified and marked according to the requirements of SP Method 2369. They are designed for the proper storage of flammable liquids, aerosol dispensers and aerosol cans in retail outlets in accordance with applicable national regulations.

## 2. TRANSPORT

### ATTENTION:

Transport the cabinet in an upright position on a pallet truck, tied and secured against slipping, until the final place of installation is reached. The transport locks in the door joints may only be removed directly at the place of installation! Inappropriate transport can lead to concealed damage to the fire protection insulation! We can only guarantee the necessary quality if the cabinet is transported to the place of its use by our specially trained staff.

### 2.1. TILTING THE CABINET

### ATTENTION:


Tilting the cabinet may only be done without jolts!

### 2.2. IN-PLANT TRANSPORT

- is also possible without transport locks (inserted as standard in the door joints), but may only be used without cabinet contents


### ATTENTION:

The model with **550 mm cabinet width** (figure 2) can only be driven in with one tine. **Pay attention to the danger of tipping!**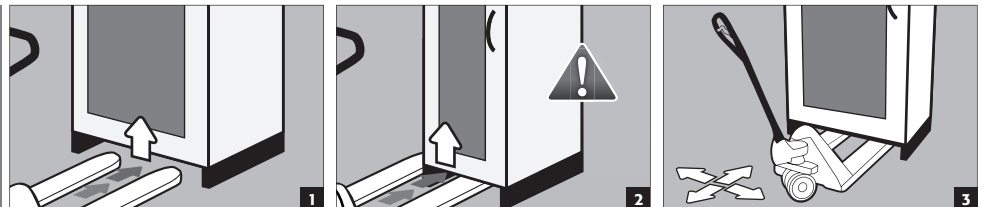


## 3. ERECTION - COMMISSIONING - FUNCTION

### 3.1. ALIGNMENT OF THE CABINETS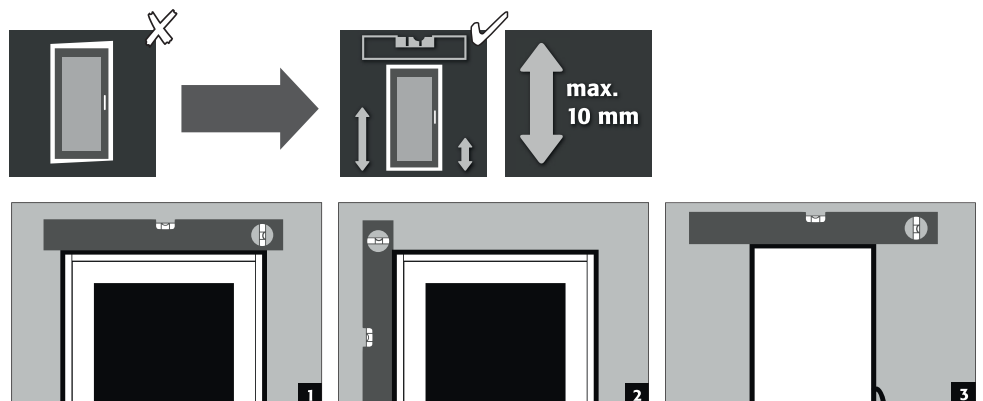

**ATTENTION:**

Door elements must not scrape against the fire prevention seals in the fold of the door when opening and closing! Doors with an automatic closing mechanism must close automatically from every position!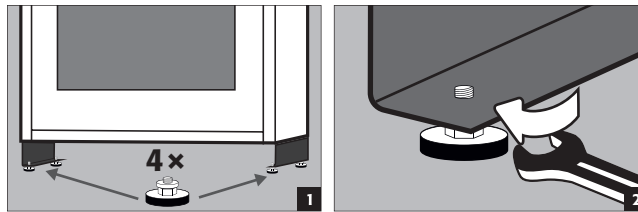

### 3.2. COMMISSIONING

Before putting into operation for the first time, the user must carry out an examination of the safety storage cabinet for possible damage, such as defective or loose sealing elements, correct alignment and perfect functioning of the door elements

**ATTENTION:**


Use the cabinet and accessories only if they are in an orderly condition.

### 3.3. INTERIOR LIGHTING (OPTIONAL)

The safety cabinets can optionally be equipped with LED lighting for the interior. For this purpose, a power supply unit for the 230 V mains connection is located on the head part of the cabinet

**Connection to the power supply**

**Equipped with:** 5 m cable including safety plug

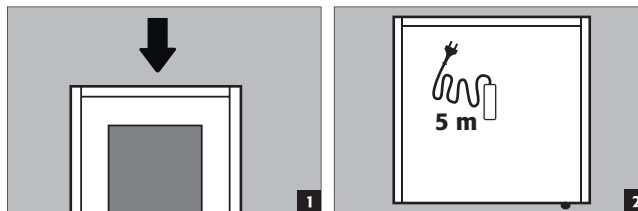**ATTENTION:**

Only use the enhanced power supply unit!

## 4. CLOSING MECHANISM

The door is permanently self-closing.


## 5. INTERIOR EQUIPMENT

### ATTENTION:


Loose parts (e.g. insertable sumps) must always be fully installed/inserted in order to ensure secure closing of the cabinet doors in case of fire.

### 5.1. BOTTOM COLLECTING SUMP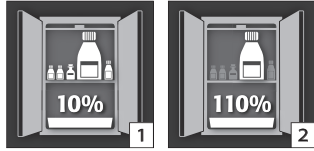

**According to SP Method 2369:** The bottom collecting sump must be able to hold a minimum volume of 10 % of all the containers stored in the cabinet **1** or at least 110 % of the volume of the largest single container **2**, whichever of these volumes is larger.

#### Use as a storage level

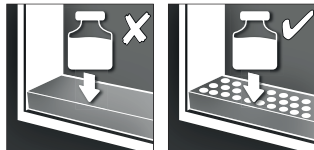
### ATTENTION


Use as a storage surface is possible only in conjunction with a perforated metal plate insert

#### Leaks:


Liquid in the sump is to be collected using suitable means. The choice of means is your own responsibility.

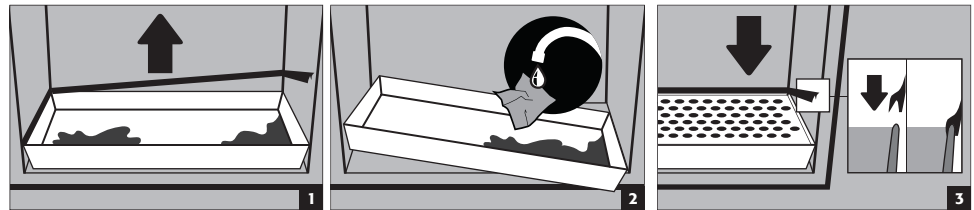### 5.2. SHELVES (HEIGHT-ADJUSTABLE)
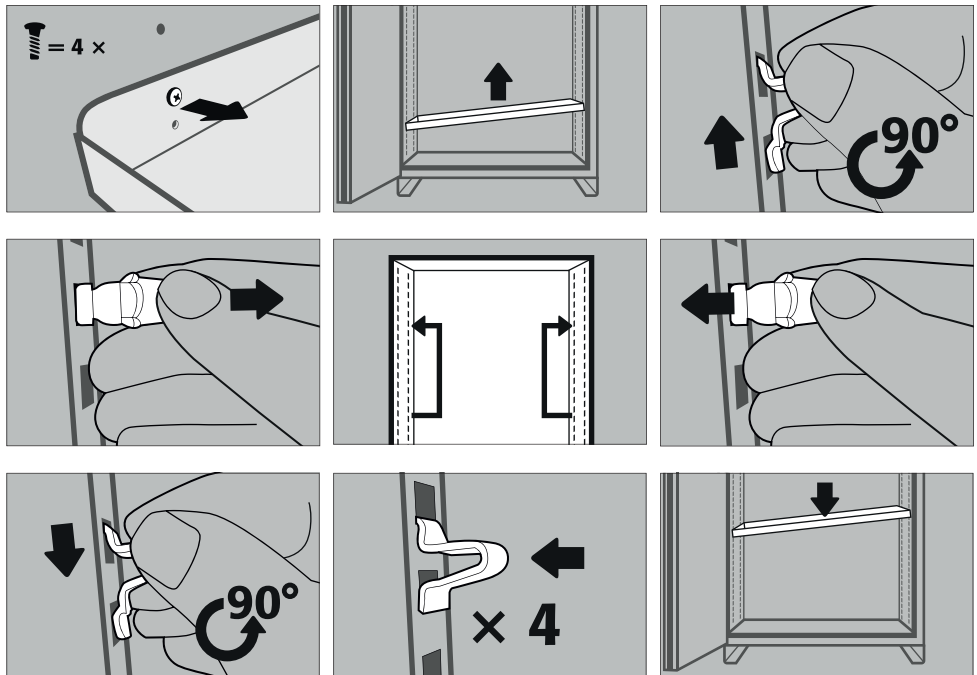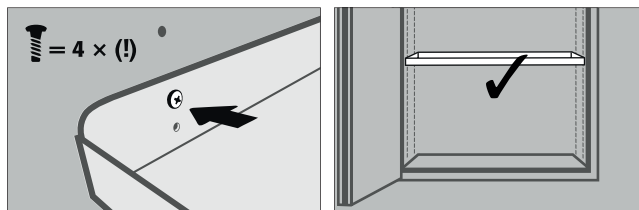
## 6. STORAGE

- Containers with aggressive chemicals (acids and alkalis) are to be kept in special cabinets for acids and alkalis or on shelves/in compartments for acids and alkalis as well as in safety storage cabinets with metal-free interior equipment.

### Cabinets prepared for connection to exhaust air system (Models EA)

- The storage of corrosive liquids can affect the function of the shut-off equipment for the air supply and exhaust.

#### Valid for all models:

Depending on the type class of the cabinet, please note the corresponding valid national regulations!

## 7. VENTILATION

### Valid for the following models

**SP30.145.085.EA.WDFW(R)**

**SP.30.195.055.EA.WDFW(R)**

**SP.30.195.085.EA.WDFW(R)**

- All models can be connected to a technical exhaust system via the exhaust connector on top of the cabinet (DN 75). Please see the respectively applicable national regulations for the specified air recirculation rate, maximum pressure losses and exhaust air ducts

#### IN GENERAL:

In general these models can be operated without technical ventilation.

## 8. FIRE - DISPOSAL

### 8.1. OPENING THE CABINET AFTER THE FIRE

#### \*ATTENTION:

**Only open the cabinet after it has cooled down. This is 6 times the fire duration!**

Depending upon the duration of the fire an ignitable vapour/air mixture may have formed; therefore remove all sources of ignition within a 10 metre radius of the cabinet before opening it

- use only spark-free tools
- open the cabinets with extreme caution

## 8.2. DISPOSAL

The models can be disposed of once they have been dismantled and the materials sorted. They are free from materials that have to be disposed of as special waste.

## 9. SAFETY CHECKS

### 9.1. GENERAL INFORMATION

As safety equipment the cabinets have to be checked for safety at least once per year. The next checking date can be taken from the check sticker on the outside of the door. This annual check can be carried out with the necessary care, and for securing your warranty claims in the case of fire, only by an authorised asecos employee (refer also to our service brochure regarding this)

**In addition, we recommend that you carry out a daily and monthly functional check:**

#### **Daily functional check**

- applies to the bottom collecting sump (collect and remove any leakages)

#### **Monthly functional check**

perfect function of the doors:

- hinges
- door closer
- correct seating and condition of the fire protection seals

In case of damage please contact your dealer in order to have the cabinet repaired using original spare parts. The cabinets can be cleaned with a mild household cleaner and a soft cloth

### 9.2. CONTACT

#### **CONTACT**


In the case of defects or complaints about our products (within and also after the warranty period), and for requesting safety checks or taking out a service contract, please contact our service hotline on:  
Tel: +49 1805 92 20 92 [info@asecos.com](mailto:info@asecos.com)

## 10. TECHNICAL DATA


		<b>SP30.145.085.EA.WDFW(R)</b> <b>SP30.145.085.NE.WDFW(R)</b>	<b>SP30.195.055.EA.WDFW(R)</b> <b>SP30.195.055.NE.WDFW(R)</b>	<b>SP30.195.085.EA.WDFW(R)</b> <b>SP30.195.085.NE.WDFW(R)</b>
Fire resistance	minutes	30	30	30
Dimensions B x T x H exterior	mm	846 × 558 × 1451	546 × 558 × 1951	846 × 558 × 1951
Dimensions B x T x H interior	mm	750 × 505 × 1285	450 × 505 × 1785	750 × 505 × 1785
Weight of empty cabinet	kg	219	230	284
Distributed load	kg/m <sup>2</sup>	621	976	671
Exhaust Air (models EA)	DN	75	75	75
Entry width of transport base	mm	683	383	683
Entry height of transport base	mm	95	95	95
Max. load of shelves	kg	75	75	75
Capacity bottom collecting sump	litres	27.5	16	27.5
Storage capacity (nominal)	litres	165	160	275
Storage capacity of 1-L-containers	pcs	153	135	255
Storage levels	pcs	3	5	5

## 11. TECHNICAL DRAWING


**SP30.145.085.EA.WDFW(R)**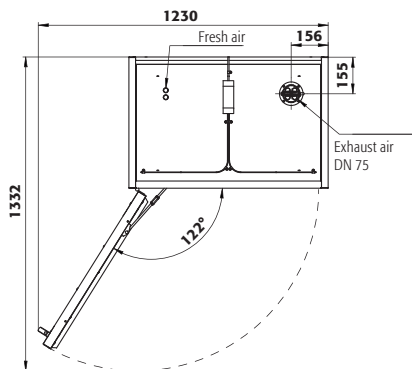


**SP30.195.055.EA.WDFW(R)**

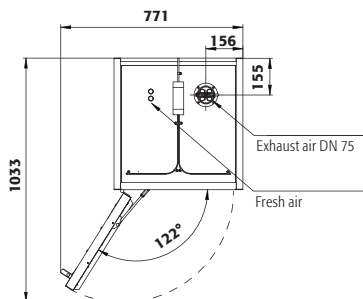
**SP30.195.085.EA.WDFW(R)**

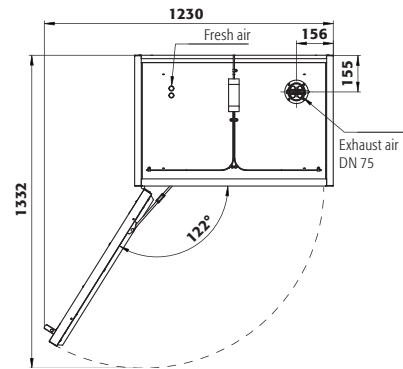
**SP30.145.085.NE.WDFW(R)**

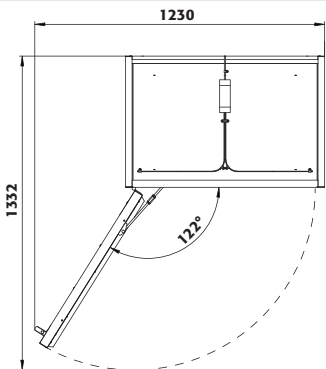
**SP30.195.055.NE.WDFW(R)**

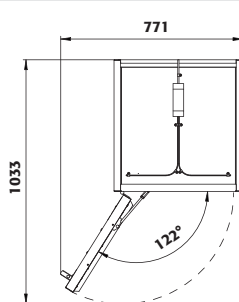
**SP30.195.085.NE.WDFW(R)**

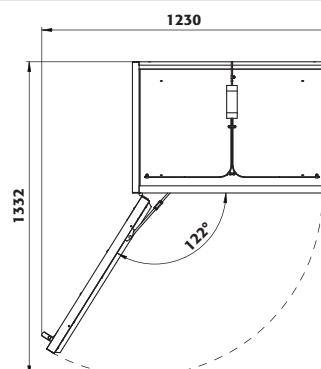
SP30.145.085.EA.WDFW(R) / SP30.145.085.NE.WDFW(R)

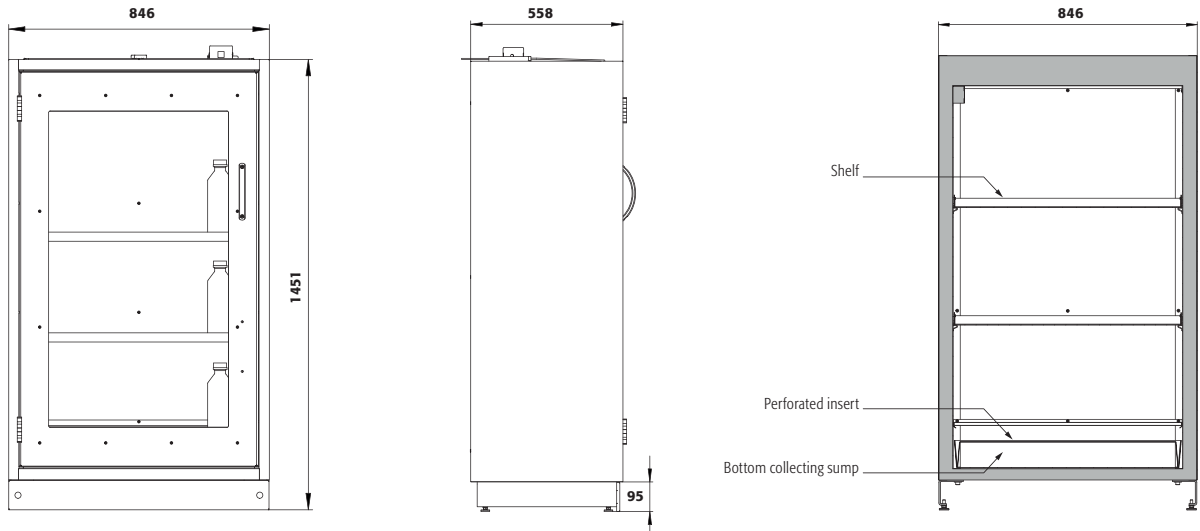SP30.195.055.EA.WDFW(R) / SP30.195.055.NE.WDFW(R)

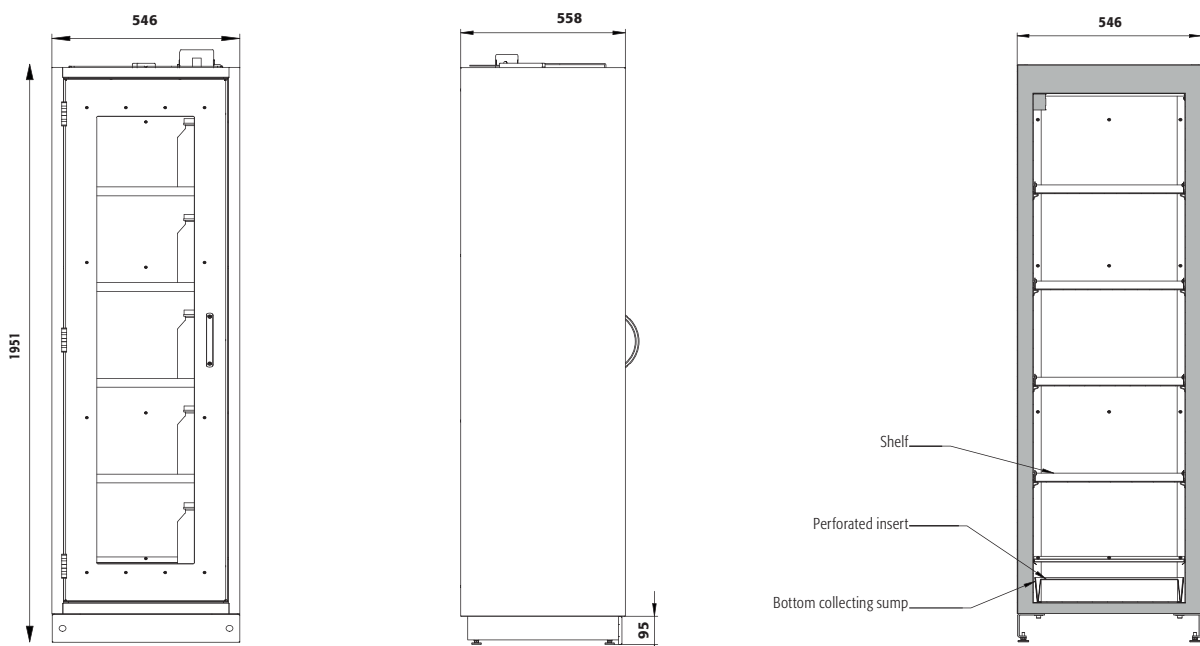SP30.195.085.EA.WDFW(R) / SP30.195.085.NE.WDFW(R)
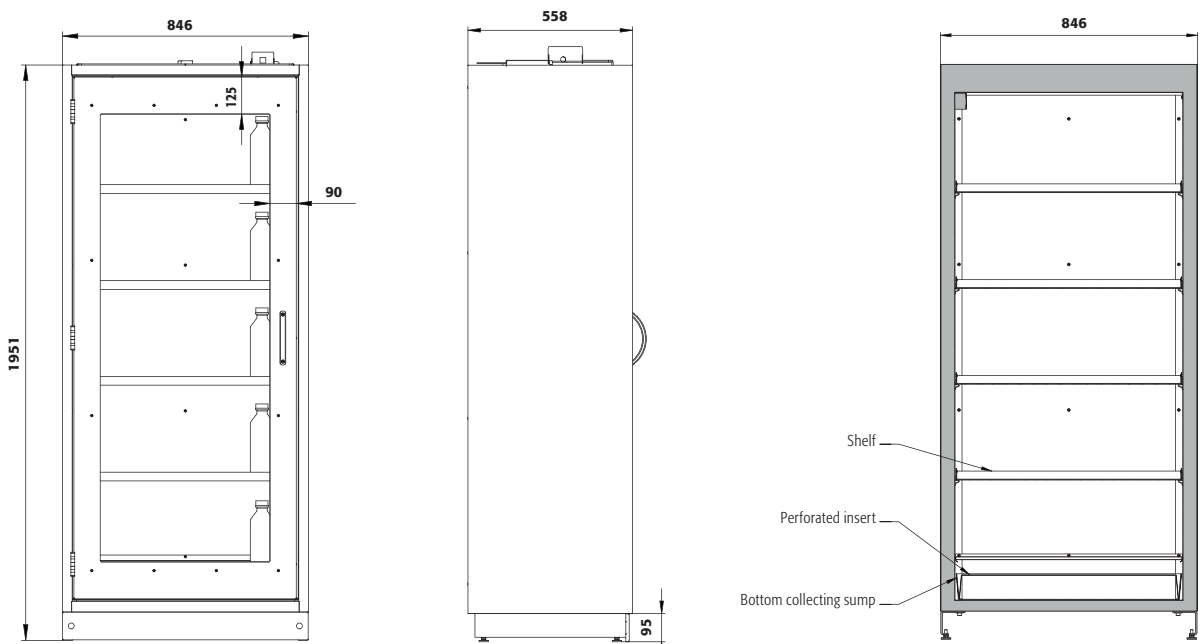
## 12. CABINET MARKING

Each cabinet model has the following three stickers on the door:

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Safety and Environmental Protection Inc.
**asecos Schweiz AG**